

Reichert SPR Concentration Analysis

Determining the Concentration of an Antigen Sample Using the SR7500DC Dual Channel Surface Plasmon Resonance (SPR) System

The concentration of an analyte can be determined directly using SPR. In this example, anti-human serum albumin mouse monoclonal (Anti-HSA) was immobilized directly to the sensor surface at a high level (greater than 1500 micro RIU). A range of antigen concentrations (human serum albumin or HSA) was injected over the surface and was expected to encompass the concentration in the unknown samples. The high surface concentration of antibody created a diffusion controlled binding response where the initial response was proportional to the concentration of the injected antigen analyte. In this example, the response at 90 seconds was plotted against the concentration of injected HSA creating a standard curve so that the concentration of HSA in unknown samples could be calculated. Known control samples were also injected.

Experimental

The experimental conditions are summarized in the following table:

- **Ligand:** Anti-HSA
- **Analyte:** HSA
- **Analyte Concentrations:** 40, 20, 10, 5, 2.5, 1.25, 0.626, 0.3125, 0.1562, 0.781, 0.0391, 0.0195nM
- **Association Time:** 3 min
- **Dissociation Time:** 3 min
- **Regeneration:** 10 mM Glycine pH 2 with 10% Glycerol

Results

The SR7500DC SPR system monitors this antibody-antigen interaction in real-time with simultaneous measurements of sample and reference channels.

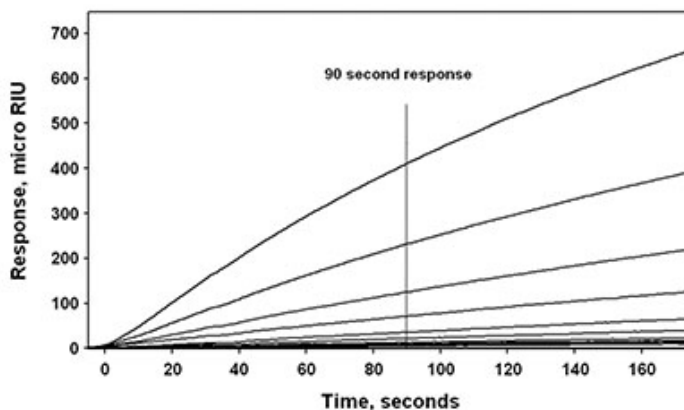


FIGURE 1

Presents normalized reference subtracted data from a series of 12 HSA injections at concentrations ranging from 0.02 to 40 nM. Between each injection the surface was regenerated with 10 mM glycine with 10% glycerol. Illustrated here is normalized response versus time plots of HSA binding to surface immobilized Anti-HSA.

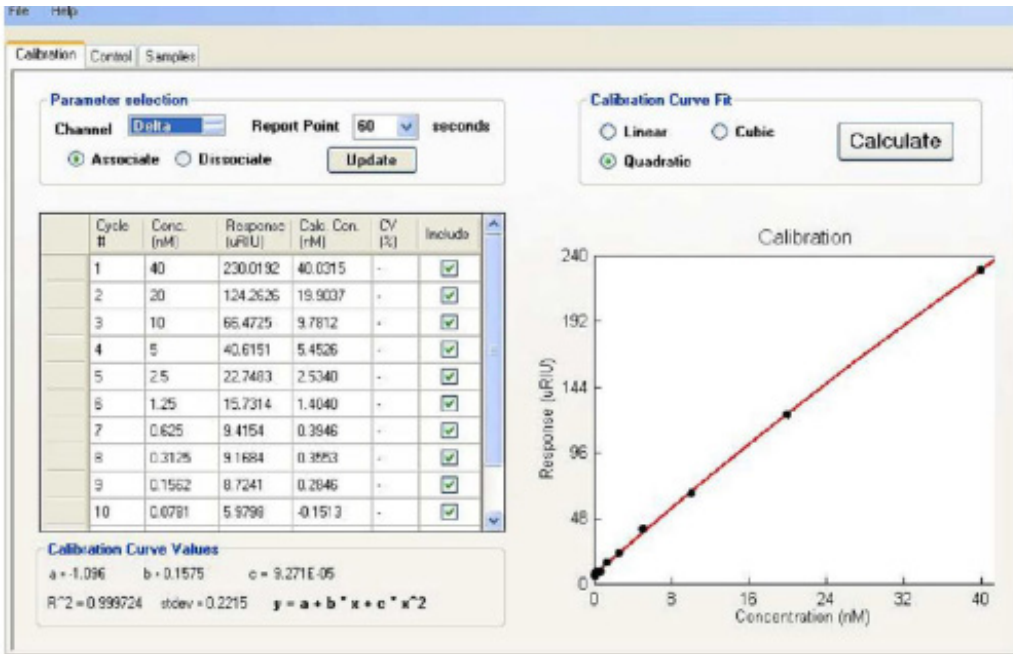


FIGURE 2

Shows the instrument response at 90 seconds for HSA standard injections versus concentration of HSA which is best fit to a quadratic. This plot can be used as a calibration curve to determine the active concentration of unknown samples. The SPR assay is a rapid direct measurement, independent of turbidity or color and does not require the use of labels or tags.

Learn more about how Reichert pushes the limits of detection and sensitivity in label-free interaction. To improve the quality of your protein interaction research, visit www.ReichertSPR.com

®/™Trademark of Reichert Technologies Company Notice: No freedom from infringement of any patent owned by Reichert or others is to be inferred. Because use conditions and applicable laws may differ from one location to another and may change with time, Customer is responsible for determining whether products and the information in this document are appropriate for Customer's use and for ensuring that Customer's workplace and disposal practices are in compliance with applicable laws and other government enactments. The product shown in this literature may not be available for sale and/or available in all geographies where Reichert is represented. The claims made may not have been approved for use in all countries. Reichert assumes no obligation or liability for the information in this document. References to "Reichert" or the "Company" mean the Reichert legal entity selling the products to Customer unless otherwise expressly noted. NO WARRANTIES ARE GIVEN; ALL IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE ARE EXPRESSLY EXCLUDED.

Corporate Office
 Reichert Technologies
 3362 Walden Avenue
 Buffalo, New York 14043, USA
 Tel: +1 716-686-4500
 Fax: +1 716-686-4555
 Toll Free USA: 1-888-849-8955
 reichertspr.lifesciences@ametek.com

European Service Center
 Carl-von-Linde Str. 42-85716
 Unterschleissheim/Munich, Germany
 Tel: +49 89 315 8911 0
 Fax: +49 89 315 891

www.reichertspr.com