


REICHERT SPR SYSTEMS

APPLICATION NOTE 6

Reichert SPR High Throughput Antibody-Antigen Analysis

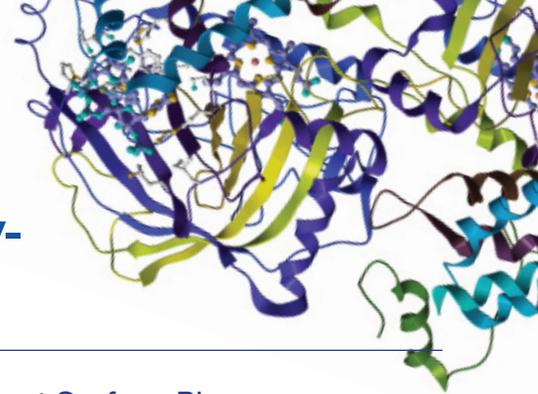

High Throughput Analysis of an Antibody-Antigen System with Reichert Surface Plasmon Resonance (SPR)

The ability to analyze a large number of samples is becoming increasingly important to researchers in academia and industry. Reichert SPR systems utilize a modular, component-based design that provides a high level of user flexibility. Part of this flexibility includes the unique ability to choose an appropriate injector that will suit the application of interest. For higher throughput applications, Reichert offers an autosampler to completely automate the injection process. The Reichert robotic autosampler can inject samples from two 96 or 384 well-plates or two 48 vial trays. Thus, the autosampler can be programmed to automatically inject samples while the researcher does other work or leaves the lab. This application note presents an antibody-antigen assay, namely bovine serum albumin (BSA) binding to anti-BSA, that includes injections from 96 different samples during a single overnight run.

Experimental

The experimental conditions are summarized in the following table:

Ligand: BSA

Analyte: Anti-BSA

Analyte Concentrations: 10, 5 and 2.5 nM

Association Time: 3 min

Dissociation Time: 4 min

Regeneration Solution: 20 mM HCL

Results

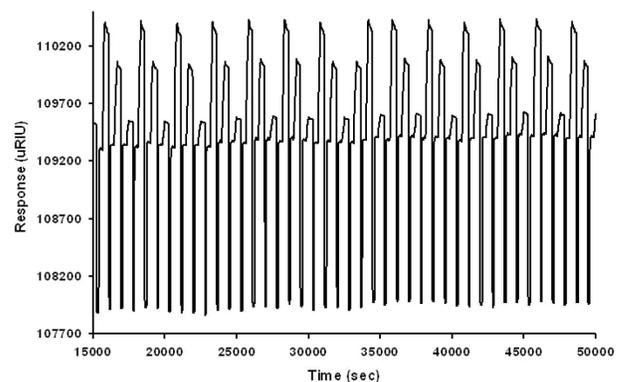
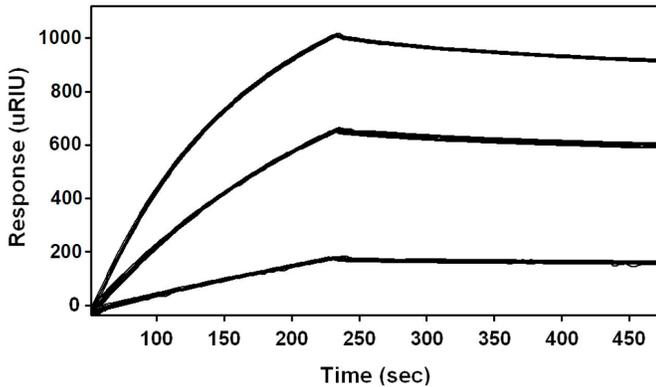


FIGURE 1

Presents a portion of the sensorgrams resulting from the 96-sample anti-BSA/BSA analysis that was programmed for a single overnight run with the Reichert robotic autosampler. Anti-BSA is injected over the sensor surface with immobilized BSA at 3 different concentrations and repeated multiple times for a total of 96 separate injections. The surface is regenerated with 20 mM HCL between each anti-BSA injection. The results indicate that the Reichert robotic autosampler injected all 96 samples without any mistakes.

**FIGURE 2**

Presents 15 overlays of each anti-BSA concentration that was injected over the sensor surface. The results show that the system is very reproducible with minimal injection-to-injection variability. This example shows that the Reichert SPR system can be used to analyze a large number of samples in a single programmed run with highly reproducible results.

Learn more about how Reichert pushes the limits of detection and sensitivity in label-free interaction. To improve the quality of your protein interaction research, visit www.ReichertSPR.com

®Trademark of Reichert Technologies Company Notice: No freedom from infringement of any patent owned by Reichert or others is to be inferred. Because use conditions and applicable laws may differ from one location to another and may change with time, Customer is responsible for determining whether products and the information in this document are appropriate for Customer's use and for ensuring that Customer's workplace and disposal practices are in compliance with applicable laws and other government enactments. The product shown in this literature may not be available for sale and/or available in all geographies where Reichert is represented. The claims made may not have been approved for use in all countries. Reichert assumes no obligation or liability for the information in this document. References to "Reichert" or the "Company" mean the Reichert legal entity selling the products to Customer unless otherwise expressly noted. NO WARRANTIES ARE GIVEN; ALL IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE ARE EXPRESSLY EXCLUDED.

Corporate Office
 Reichert Technologies
 3362 Walden Avenue
 Buffalo, New York 14043, USA
 Tel: +1 716-686-4500
 Fax: +1 716-686-4555
 Toll Free USA: 1-888-849-8955
reichertspr.lifesciences@ametec.com

European Service Center
 Carl-von-Linde Str. 42-85716
 Unterschleissheim/Munich, Germany
 Tel: +49 89 315 8911 0
 Fax: +49 89 315 891

www.reichertspr.com